


1,3-[Bis(diphenylphosphino)propane] nickel(II) chloride

Product number: 416 | Alternative name: NiCl₂(dppp) | CAS number: 15629-92-2 | C₂₇H₂₆Cl₂NiP₂ | Molecular weight: 542.04

NiCl₂(dppp) is an air stable, dark red or purple square planar complex. NiCl₂(dppp) is a very useful catalyst in preparation of C-C bonds via coupling with Grignard (Kumada reaction) and zinc (Negishi reaction) reagents.

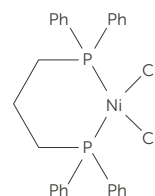


NICKEL-CATALYZED CROSS-COUPLING WITH ZINC AND GRIGNARD REAGENTS¹⁻¹²




R, R' = alkyl, alkenyl, aryl heterocyclic
 X = halides, dialkylcarbonates, arenesulfonates, heteraromatic ethers, phosphates
 Y = Zn, Mg

PREPARATION OF HIGHLY CONDUCTIVE, REGIOREGULAR POLYMERS^{15, 16}

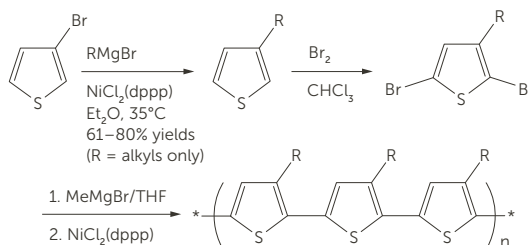


R = alkyls, alkoxys, esters HT regioregularity 91 to >99% Mw range 12,000 up to 50,000 PDIs = 1.1 to 1.78

FACILE AND SELECTIVE DEALLYLATION OF ETHERS, AMINES, AND AMIDES^{13,14}

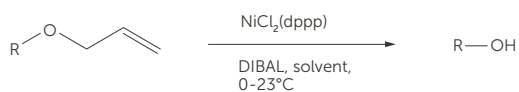


R = alkyl, alkenyl, aryl, carbohydrates 71 - 97% yield

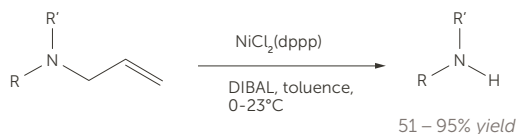

R = H, alkyl
 R' = H, Me, Bn, CH₂CH=C(Me)₂, allyl, p-tolSO₂, C(O)R'
 R/R' = heterocycles

HYDROBORONYLATION OF THIOALKYNES¹⁷

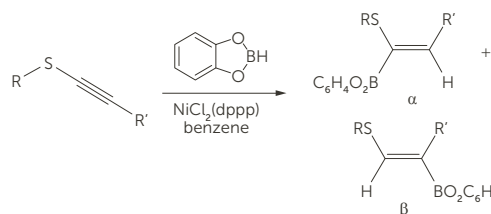

a: R = Et, R' = Me 93 - 97% yields
 b: R = Ph, R' = H β/α: 98:2

References

- ¹ Erdelmeier, I. and Gais, H.-J., *J. Am. Chem. Soc.*, **1989**, *111*, 1125.
- ² Tagat, J.R., McCombie, S.W., Barton, B.E., Jackson, J., and Shortall, J., *Bioorg. Med. Chem. Lett.*, **1995**, *5*(18), 2143.
- ³ Tamao, K., Kiso, Y., Sumitani, K., and Kumada, M., *J. Am. Chem. Soc.*, **1972**, *94*, 9268.
- ⁴ Sengupta, S., Leite, M., Raslan, D.S., Quesnelle, C., and Snieckus, V., *J. Org. Chem.*, **1992**, *57*, 4066.
- ⁵ Organ, M.G. and Murray, A.P., *J. Org. Chem.*, **1997**, *62*, 1523.
- ⁶ Hopf, H. and Theurig, M., *Angew. Chem. Chem. Int. Ed.*, **1994**, *33*(10), 1099.
- ⁷ Bigas, A.F. and Johnston, R.A.W., *J. Chem. Soc., Perkins Trans. 1*, **2000**, 1735.
- ⁸ Poree, J.-H., Clavel, A., Betzer, J.-F., Pancrazi, A., and Ardisson, J., *Tetrahedron Lett.*, **2003**, *44*, 7553.
- ⁹ Kitamura, M. and Narasaka, K., *Org. Syn.*, **2001**, *78*, 42.
- ¹⁰ Cho, C.-H., Sun, M., Seo, Y.-S., Kim, C.-B., and Park, K., *J. Org. Chem.*, **2005**, *70*, 1482.
- ¹¹ Sofia, A., Karlstrom, E., Itami, K., and Backvall, J.-E., *J. Org. Chem.*, **1999**, *64*, 1745.
- ¹² Mans, D.M. and Pearson, W.H., *J. Org. Chem.*, **2004**, *69*, 6419.
- ¹³ Taniguchi, T. and Ogasawara, K., *Angew. Chem. Int. Ed.*, **1998**, *37*(8), 1136.
- ¹⁴ Taniguchi, T. and Ogasawara, K., *Tetrahedron Lett.*, **1998**, *39*, 4679.
- ¹⁵ McCullough, R.D. and Lowe, R.D., *J. Chem. Soc., Chem. Commun.*, **1992**, 70.
- ¹⁶ Amarasekara, A.S. and Pomerantz, M., *Synlett*, **2003**, 2255.
- ¹⁷ Gridnev, I.D., Miyaura, N., and Suzuki, A., *Organometallics*, **1993**, *12*, 589.

FOR MORE INFORMATION

Please call your Regional Customer Service Center today to learn what Entegris can do for you. Visit entegris.com and select the [Contact Us](#) link to find the customer service center nearest you.

TERMS AND CONDITIONS OF SALE

All purchases are subject to Entegris' Terms and Conditions of Sale. To view and print this information, visit entegris.com and select the [Terms & Conditions](#) link in the footer.



Corporate Headquarters

129 Concord Road
Billerica, MA 01821
USA

Customer Service

Tel +1 952 556 4181
Fax +1 952 556 8022
Toll Free 800 394 4083

Entegris®, the Entegris Rings Design®, and other product names are trademarks of Entegris, Inc. as listed on entegris.com/trademarks. All third-party product names, logos, and company names are trademarks or registered trademarks of their respective owners. Use of them does not imply any affiliation, sponsorship, or endorsement by the trademark owner.

©2018-2020 Entegris, Inc. | All rights reserved. | Printed in the USA | 9000-11206TAN-1120