

AccuSizer A2000 MPA Microplate Particle Analyzer

*Integrated microplate automated
particle analyzer*

The AccuSizer A2000 MPA is an automated microplate reader for performing particle count and size analysis. One or two 96 well microplates are loaded into the system and the AccuSizer single particle optical sizing (SPOS) technique generates results in particles/mL from 0.5 – 400 µm. Automated USP 787 calculations are reported along with detailed size distribution, if desired.

The system (images at right) consists of the following components:

- LE400-05 light extinction/scattering sensor
- Size range 0.5 – 400 µm
- Pulse height analyzer/counter
- Autosampler to transfer sample from microplate wells through sensor

Measurement procedure:

- Microplate is loaded onto autosampler
- User sets up measurement protocol and plate configuration
- Analysis started
- Sample needle moves to first position
- Sample mixed by push/pull, if desired
- Small volume aspirated through sensor for analysis
- Sample returned to well
- Needle moves to two washing stations to clean outside and inside
- Report(s) generated
- Repeat

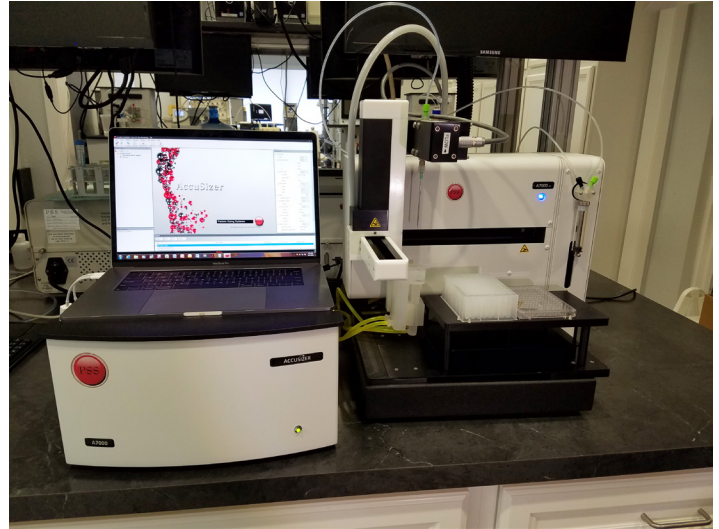


Figure 1

USP 788 - 04/27/2018 11:04:19 – Page 1 of 1

Entegris

Operator: admin ad admin
 Sample Volume: 0.25 mL
 Flow rate: 30 mL
 Report Date/Time: 04/27/2018, 11:04

Standard: USP 788 2014
 Protocol: 1 um Standard
 Project: 1 um Standard

Sample: USP 787
 Comment:

Sample	Run Date/Time	Sample Volume (mL)	≤10 µm (#/mL)	>25µm (#/mL)
Tube 1 Rep. 2	10/04/2016 22:55	0.25 mL	1914	1704

TEST Criteria	RESULT
#mL > 10 µm AND #mL > 25 µm (PASS)	FAIL

FEATURES

- Mount one or two 96 well plates
- Well volume = 1 mL
- Minimum sample volume = 180 µL
- Analysis time including cleanup = 2 minutes
- Automated USP 787 calculations
- Data exported to Excel or LIMS
- Sample conservation

FOR MORE INFORMATION

Please call your Regional Customer Service Center today to learn what Entegris can do for you. Visit entegris.com and select the [Contact Us](#) link to find the customer service center nearest you.

TERMS AND CONDITIONS OF SALE

All purchases are subject to Entegris' Terms and Conditions of Sale. To view and print this information, visit entegris.com and select the [Terms & Conditions](#) link in the footer.



Corporate Headquarters

129 Concord Road
Billerica, MA 01821
USA

Customer Service

Tel +1 952 556 4181
Fax +1 952 556 8022
Toll Free 800 394 4083

Entegris®, the Entegris Rings Design®, and other product names are trademarks of Entegris, Inc. as listed on entegris.com/trademarks. All third-party product names, logos, and company names are trademarks or registered trademarks of their respective owners. Use of them does not imply any affiliation, sponsorship, or endorsement by the trademark owner.

©2018-2019 Entegris, Inc. | All rights reserved. | Printed in the USA | 7127-10464TAN-0419